



Pulse EPR console for spectroscopy and imaging



Connecting Spectrometers to People

Boris Epel, Igor Gromov, Stefan Stoll

Visit www.specman4EPR.com for recent updates and developments. SpecMan4EPR is distributed by Scientific Software Services. Contact Reef Morse at reefmorse@scientific-software.com for information regarding purchase options and support.

Every pulse spectrometer has a console for pulse generation and detection of time domain signals. We present an affordable pulse EPR console built from general purpose components that works as a single unit. Major components of the console are:

- ◆ **SpecMan4EPR** – versatile control and acquisition software for pulse EPR experiments¹.
- ◆ **PulseBlasterESR-Pro 400 MHz**, SpinCore Technologies, Inc (www.spincore.com) – universal pulse programmer with 2.5 ns time resolution.
- ◆ **AP235 averager**, Agilent Technologies (www.acqiris.com) – 0.5 GS/s (2 ns resolution) for 2 channel acquisition with 133MB/s data transfer rate.
- ◆ National Instruments **DAQmx**, **VISA** and **GPIB** devices; **LPT**, **COM** ports; TCP-IP interfaced third-party devices.

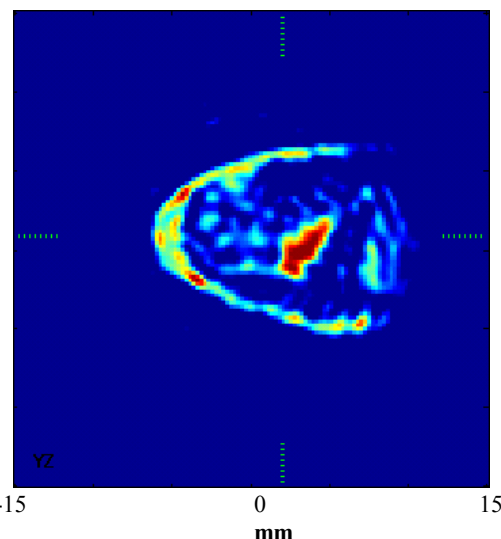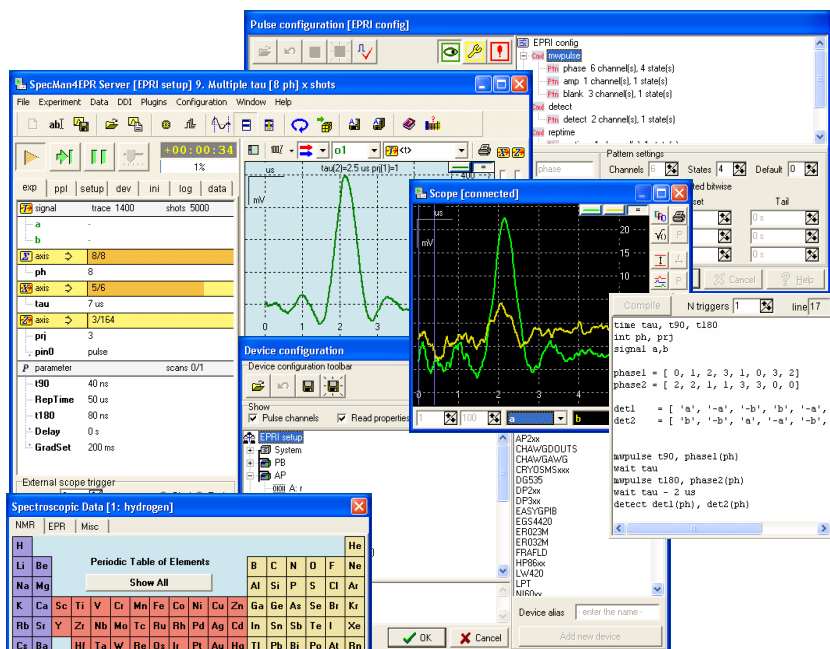

Slice of a 3D EPR image (300 μm resolution) of tumor bearing mouse leg obtained using SpecMan4EPR. 250 MHz pulse imager; multi-B ESE algorithm; 5 G/cm gradient; ~ 1200 projections; partially deuterated OX063 spin probe; 50 min acquisition time. Data are a courtesy of H.J.Halpern group, University of Chicago.

Front end windows and dialogs of the SpecMan4EPR

- ◆ Complete pulse console inside one computer case. Single system for all kinds of spectrometers.
- ◆ Device-independent pulse programming language; acquisition of multiple time traces during single pulse sequence; minimal reprogramming time.
- ◆ Four-dimensional experiments; linear, logarithmic or table-based definition of ANY device or experiment parameter.
- ◆ Remote control over LAN or Internet; **TCP-IP interface to LabView™ modules.**
- ◆ In-scope Fourier transformation and baseline correction; time-trace baseline subtraction.
- ◆ **NEW!** Queue your experiments in Batch mode; sequential or time-based execution of the queue.
- ◆ **NEW!** 0.4 GB data arrays for high resolution imaging.
- ◆ **NEW!** User-defined information fields store important experimental parameters with your data.

¹Epel B. *et al.*, J. Magn. Reson., **164**, 78 (2003); Epel B. *et al.*, Concepts in Magnetic Resonance, **26B**, 36 (2005)